

Dynamic Gas-Solid Surface Interactions for Dry Reforming of Methane via In-Situ DRIFTS Analysis

Md Monir Hossain, Ruigang Wang (rwang@msu.edu), Department of Chemical Engineering and Materials Science, Michigan State University, East Lansing, MI 48824, United States.

In this note: monitoring surface species on a nanorod catalyst for dry reforming of methane.

Dry reforming of methane (DRM) is a key process for converting CH_4 and CO_2 into valuable syngas (CO and H_2) [1–3]. Understanding gas-solid surface interactions during DRM is essential for the rational design of catalysts.

Here, in-situ Diffuse Reflectance Infrared Fourier Transform Spectroscopy (DRIFTS) using the Harrick Praying Mantis™ accessory and a High-Temperature Reaction Chamber (HVC-DRM-5) is employed to track surface species on a 1.5 wt% Ni–0.5 wt% Ru/ CeO_2 nanorod (NR) catalyst over a temperature range of 50 °C to 450 °C.

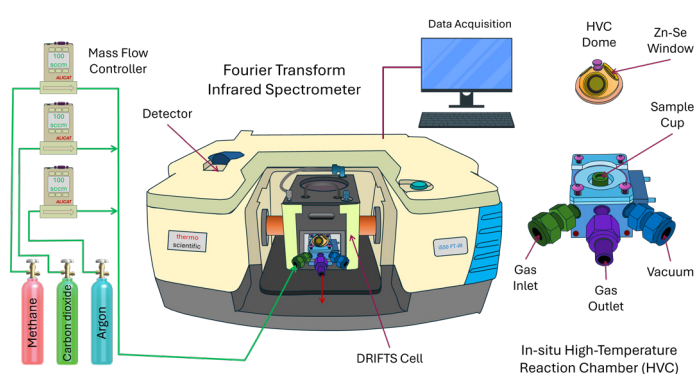


Figure 1: Experimental setup for In-Situ DRIFTS using a Harrick Praying Mantis Cell and High-Temperature Reaction Chamber.

Experimental

Catalyst Preparation: The 1.5 wt% Ni–0.5 wt% Ru/ CeO_2 NR catalyst was synthesized via a one-pot hydrothermal method, ensuring uniform dispersion of Ni–Ru on the CeO_2 NR supports.

Instrumentation: In-situ DRIFTS was performed with the Harrick Praying Mantis diffuse reflection accessory along with a High-Temperature Reaction Chamber, integrated into a Thermo Fisher Nicolet iS50 FTIR system (Fig. 1), providing precise temperature and gas flow control for real-time monitoring of surface gas-solid reactions.

Reaction Conditions: A CH_4/CO_2 gas mixture (29:71 $\approx$ 0.4 ratio) was fed at a total flow rate of 100 sccm. The temperature was ramped from 50 °C to 450 °C at 10 °C/min, with spectra collected every 50 °C after stabilizing the gas flow and temperature for 15 min.

Results and Discussion

The DRIFTS spectra of the 1.5 wt% Ni–0.5 wt% Ru/ CeO_2 NR catalyst during DRM reveal several changes in surface reaction intermediates (Fig. 2). The key observations are summarized below.

- At 50 °C, CO_2 adsorbs on the catalyst surface as various carbonate species (Fig. 2(a)) [4,5].
- Beginning at 300 °C, the bidentate carbonate species ($\sim 1020 \text{ cm}^{-1}$) begin to convert into monodentate carbonate species ($\sim 1078 \text{ cm}^{-1}$), indicating the progression of the reaction [5].
- In the CO region (Fig. 2(b)), several peaks emerge [6], with pronounced changes observed from 300 °C onward. As the temperature increases, the intensity of the adsorbed CO peaks grows, and gaseous CO peaks (~ 2120 and 2170 cm^{-1}) appear at 450 °C, signifying weakened CO adsorption and the release of CO gas.

These observations provide valuable insight into molecular bond transformations involved in the reaction mechanism. Further analysis of in-situ DRIFTS data will offer a deeper understanding of gas-solid interactions and gas-gas conversions at elevated temperatures.

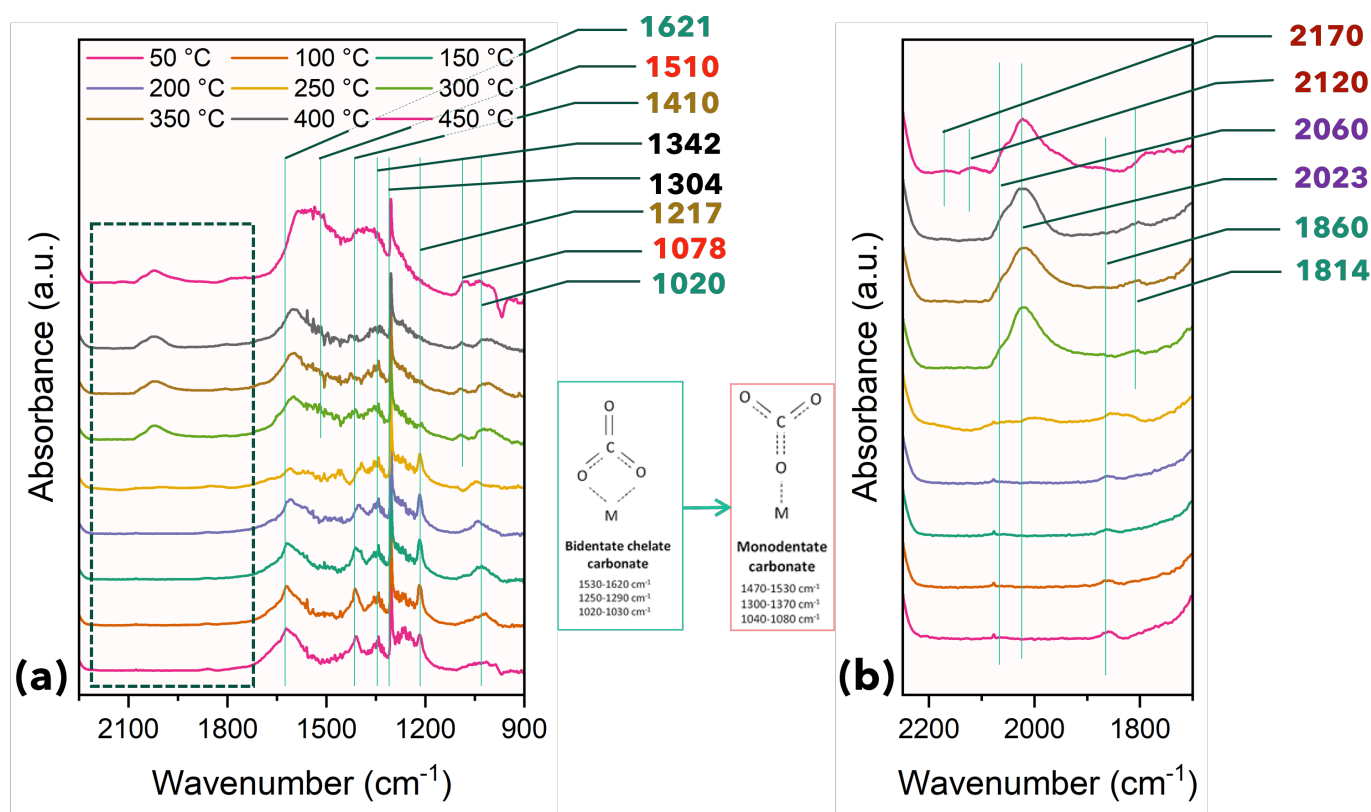


Figure 2: In-Situ DRIFTS spectra for DRM at varying temperatures ($\text{CH}_4/\text{CO}_2 = 29/71$ sccm ≈ 0.4 ratio; total flow rate: 100 sccm).

Conclusion

The in-situ DRIFTS investigation with the Harrick Praying Mantis setup revealed the evolution and transformation of surface species during DRM on Ni-Ru/CeO₂ NRs. The findings demonstrate the power of in-situ DRIFTS's ability to elucidate catalytic mechanisms and highlight the promising performance of this catalyst for DRM.

Acknowledgement

This project is supported by the U.S. National Science Foundation (2427238 and 2427213).

References

1. Ahasan, M. R.; Hossain, M. M.; Wang, R. Dielectric Barrier Discharge Reactors for Plasma-Assisted CO₂ and CH₄ Conversion: A Comprehensive Review of Reactor Design, Performance, and Future Prospects. *Energy Technology* 2024, 13 (4), 2401177.
2. Monir Hossain, M.; Robayet Ahasan, M.; Wang, R. Influence of Catalyst Shape on Plasma-Assisted Dry Reforming of Methane:

A Comparative Study of Ni-CeO₂ Nano-Cubes and Nano-Octahedra. *Chemical Engineering Journal* 2024, 496, 154193.

3. Hossain, M. M.; Ahasan, M. R.; Wang, R. Exploring the Versatile Roles of Microwave Energy in Catalyst Synthesis, Treatment, and Catalytic Performance. *Fuel* 2025, 400, 135684.
4. Vakili, R.; Gholami, R.; Stere, C. E.; Chansai, S.; Chen, H.; Holmes, S. M.; Jiao, Y.; Hardacre, C.; Fan, X. Plasma-Assisted Catalytic Dry Reforming of Methane (DRM) over Metal-Organic Frameworks (MOFs)-Based Catalysts. *Appl Catal B* 2020, 260, 118195.
5. Cárdenas-Arenas, A.; Quindimil, A.; Davó-Quiñonero, A.; Bailón-García, E.; Lozano-Castelló, D.; De-La-Torre, U.; Pereda-Ayo, B.; González-Marcos, J. A.; González-Velasco, J. R.; Bueno-López, A. Isotopic and in Situ DRIFTS Study of the CO₂ Methanation Mechanism Using Ni/CeO₂ and Ni/Al₂O₃ Catalysts. *Appl Catal B* 2020, 265, 118538.
6. Li, Y.; Qin, T.; Wei, Y.; Xiong, J.; Zhang, P.; Lai, K.; Chi, H.; Liu, X.; Chen, L.; Yu, X.; Zhao, Z.; Li, L.; Liu, J. A Single Site Ruthenium Catalyst for Robust Soot Oxidation without Platinum or Palladium. *Nature Communications* 2023 14:1 2023, 14 (1), 1–11.

Contact us today

Reach out to us any time for more information on the application described in the text above or visit our website [specac.com](https://www.specac.com) to read more and make an enquiry.

United Kingdom
sales@specac.co.uk
+44 (0) 1689 892 902

United States
sales@specac.com
+1 800-248-3847

China
cn@specac.com

Singapore
sg@specac.co.uk

